GIMA

Codice 24035 Centrifuga XC-2000

Code 24035 XC-2000 Centrifuge

XC-2000 CENTRIFUGA

Istruzioni

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XC-2000 CENTRIFUGE

Operation

1. Open the lid. (The unit is designed to automatically stop, when the lid is open.)
2. Place supplied plastic tube holder shells into the centrifuge holes. Holders are provided in two sizes, 15ml & 10ml. Use the appropriate size holder for your tubes.
3. Insert the centrifuge tubes that you want to spin in a manner to balance out the specimen tubes' weight. In order to get a smooth running centrifuge, always insert even number of tubes on opposite sides of centrifuge with approximately the same amount of fluid in each tube.
4. Using the one touch switch panel: push the "switch" button to get to the time function. Push the + or - button to adjust the time setting: 1 minute to 60 minutes. Push the "switch" button to get to the speed function. Push the + or - button to adjust the speed setting: 1000 rpm to 4000 rpm in 500 rpm increments.
5. Push the start button to begin program. This centrifuge has a memory setting and will run the previous program.
6. If the tube is too short for the tube holder shell, you may use some of the soft foam provided by the manufacturer. Please make sure that specimen tubes are snug into holder before starting centrifuge.

Safety Instructions

- This machine is designed for indoor use on a flat surface.
- Make sure the specimen tubes are equal and balanced. Unbalanced tubes will cause the machine to shake and not provide a smooth spinning motion.
- Make sure specimen tubes are completely sealed before turning on the centrifuge.
- Allow centrifuge to come to a complete and full stop before removing the specimens.
- Disconnect from power source before attempting to clean the inside of this machine.
- Never move the centrifuge while it is running.

Specifications

Voltage: 220v-50hz

Speed: Approx. 1000-4000 rpm (increased in increments of 500 rpm's)

Time: 1-60 min.

Capacity: 15ml or 10 ml x6

Max.RCF: 1790g

Power: 80w

Maintenance

Always keep the outside and inside of the machine clean. A dustcover is recommended while it is not in use.

If you have problems operating the centrifuge, please call your dealer or authorized service provider for help. Do not attempt to fix the machine yourself.